

Thermo Scientific ISQ Series Single Quadrupole GC-MS

Advanced GC-MS systems designed for continuous high-throughput operation

With 50 years of GC-MS innovation, the Thermo Scientific™ ISQ™ Series continues to deliver quality design and leading technology in single quadrupole mass spectrometry. The series consists of the ISQ QD GC-MS and ISQ LT GC-MS models.

For budget-minded lab owners seeking a reliable replacement for older GC-MS or stand-alone GCs, the ISQ QD GC-MS models represent robust solutions with quality design, MS operational simplicity, and proven dependability.

For high-throughput, analytically demanding laboratories, requiring the utmost sensitivity and unstoppable productivity, the ISQ LT GC-MS offers future-proof investment into the leading technology, simply intelligent GC-MS platform. Unlike other systems, the ISQ LT GC-MS does not need to be vented to exchange the ion source and provides utmost sensitivity, low detection limits and unparalleled flexibility.



Thermo
SCIENTIFIC

ISQ Series GC-MS Specifications

Mode (Standard)

- Electron Impact Ionization (EI), with full scan (FS), SIM, and FS/SIM simultaneous within sample injection

Modes (Optional)

- AutoSIM and timed acquisition (t-SIM) mode (standard on ISQ LT and ISQ QD 300, optional on ISQ QD)
- Chemical ionization, including PCI, NCI, Thermo Scientific PPINICI and EI/PCI/NCI combination (ISQ LT only)

Ion Source Type

- Thermo Scientific™ ExtractaBrite™ off-axis ion source, with repeller, source lenses, RF lens and dual filaments in all ionization modes
- Programmable to 350 °C

Source Access

- Vacuum probe interlock to remove entire source or change ionization modes without venting (standard and available on ISQ LT model only)
- Combination EI/PCI/NCI source used without interchange

Mass Filter

- Dual-stage mass filter with off-axis ion guide pre-filter for noise reduction and solid, homogeneous non-coated, maintenance-free quadrupole rods
- Independently heated zones for: transfer line, ion source, and ion optics/mass analyzer
- Fast quadrupole scanning, up to 20,000 u/s

Mass Stability

- Better than 0.1 u/48 hours/ $\Delta T \leq 2$ K

Detector

- Triple off-axis DynaMax XR detection system, with off-axis 10 kV dynode, discrete dynode electron multiplier and electrometer, linear from 0 to 68 μ A
- Electrical dynamic range $> 10^9$

Mass Range

- 1.2–1100 u with unit mass resolution

Acquisition Rate

- Ability to acquire more than 240 scans/s in SIM
- Ability to acquire more than 97 scans/s in FS when scanning over a range of 125 u

Pumping Systems

- Standard capacity (66 L/s) turbomolecular pump (ISQ QD GC-MS)
- Extended capacity (300 L/s) turbomolecular pump (ISQ LT GC-MS and ISQ QD 300 GC-MS)
- Mechanical rotary vane 3.3 m³/h oil pump
- Optional oil-free scroll pump

CI Reagent Gas Capabilities

- Software-switchable dual reagent gas with digital flow control
- Ammonia, methane, isobutane and CO₂ presets

Electron Energy

- Adjustable from 0 eV to 150 eV

Emission Current

- Up to 350 μ A

Transfer Line Temperature

- Up to 400 °C

Gas Chromatograph (Thermo Scientific™ TRACE™ 1300 Series GC)

Please refer to GC and autosampler specification sheets for additional and up-to-date details

- TRACE 1300 GC: Intuitive, single-button start/stop for ease of use with minimal local instrument interaction
- TRACE 1310 GC: Complete icon-driven touch screen user interface for direct local instrument control

Autosamplers

- AI/AS 1310, TriPlus RSH, TriPlus 100 LS, TriPlus 300 HS, and more

Instant Connect Modules

- User-installable injector or detector assembly can be installed in less than 2 minutes
- 0.001–1000 kPa digitally controlled carrier gas with gas saver and septum purge
- Detector Fast Data Acquisition Rate: up to 300 Hz

Oven Performance

- Operating range: Ambient +3 °C to 450 °C
- Cryogenic Option Minimum Temperature: -100 °C with LN₂; -50 °C with liquid CO₂
- Maximum heating rate: 125 °C/min
- Oven cool-down (22 °C ambient): 450 °C to 50 °C in <4 min

GC Analytical Performance

- Retention Time Repeatability: <0.0008 min
- Peak Area Repeatability: <0.5 % RSD

Hydrogen Kit Option

- Optional Hydrogen Kit (includes hydrogen sensor and ion volume) and extended capacity turbomolecular pump required for use with and standard installation specifications performed with hydrogen carrier gas on the TRACE 1300/1310 GC with the ISQ Series MS

Instant Connect Helium Saver Module

- Compatible with any ISQ Series MS model connected to a TRACE 1300 Series GC
- Save helium during the analytical run, and when instrument is idle
- Analytical conditions remain the same; methods remain intact
- Realize significant savings in helium supply throughout the lifetime of your GC or GC-MS instrument

Direct Sample Probe System Option (ISQ LT Only)

- Switch to probe in <3 min with GC undisturbed
- Available in two styles: rapid heating filament Direct-Exposure Probe (DEP, capable of flash vaporization or pyrolysis at up to 1600 °C) or slower volatilization Direct-Insertion Probe (DIP, capable of accommodating powders and solid samples in a quartz or aluminum crucible) up to 450 °C

Microfluidics Options for TRACE 1300/1310 GC

- NoVent for easier column replacement without venting the system
- Dual Detector kit for splitting column effluent to two detectors (including MS)
- Consists of low-volume, highly inert SilFlow™ technology with finger-tight connectors

Data System Software and Options

- Thermo Scientific™ TraceFinder™ Software for routine GC-MS quantitation
- Thermo Scientific™ Xcalibur™ Data system
- Thermo Scientific™ Dionex™ Chromeleon™ 7.2 CDS for chromatographers using MS
- TraceFinder Software for Environmental and Food Safety
- TraceFinder Software for Clinical Research
- TraceFinder Software for Forensic Toxicology
- Instrument control and data connection via Ethernet
- Computer supplied with instrument equipped with three Ethernet (8P8C RJ-45) ports
- Commercial mass spectral library options, including
 - NIST and Wiley libraries
 - Mass Spectral and GC Data of Drugs, Poisons, Pesticides, Pollutants and their Metabolites (Maurer-Pfleger-Weber library)

Standard Installation Specifications* ISQ LT and ISQ QD 300 (ISQ QD)

Mode/Concentration	He	H ₂
In EI mode, 1 μ L of 1 pg/ μ L octafluoronaphthalene (OFN) will produce the following minimum signal-to-noise for <i>m/z</i> 272 when scanning 50–300 u	1,500:1 (600:1)	100:1
In PCI mode, 1 μ L of 100 pg/ μ L benzophenone will produce the following minimum signal-to-noise for <i>m/z</i> 183 when scanning 80–230 u using methane reagent gas	300:1	300:1
In NCI mode, 2 μ L of 100 fg/ μ L of OFN will produce the following minimum signal-to-noise for <i>m/z</i> 272 when scanning 50–300 u using methane reagent gas	600:1	600:1

Typical Specifications

In EI SIM mode, with helium carrier gas and the AI/AS 1310, TriPlus 100 LS, or TriPlus RSH autosampler (required and configured for liquid injections), eight sequential 100 fg OFN splitless injections monitored for *m/z* 272 produce the following instrument detection limit (IDL), calculated from the chromatographic peak area with 99% confidence interval: **IDL \leq 10 fg (ISQ LT GC-MS and ISQ QD 300 GC-MS); IDL \leq 24 fg (ISQ QD GC-MS)**

System Dimensions/Weights

Total width of the connected GC/MS system is 80 cm (31 in). Allow 16 cm (6 in) of clearance behind the instrument (32 cm if using autosampler). Additional space should be allotted for data system and printer.

Mass Spectrometer

(height \times width \times depth)

44 \times 33 \times 63 cm (17.5 \times 13 \times 24.5 in)

Weight: 43 kg (94 lbs)

TRACE 1300/1310 GC

45 \times 44 \times 60/67 cm (18 \times 17 \times 24/26 in)

Weight: 35 kg (77 lbs)

* He (H₂) standard specifications are performed using a 15 (30) m \times 0.25 mm ID \times 0.25 μ m System Qualification Column (SQC). The installation specifications are performed with either He or H₂ but not both.

www.thermoscientific.com

©2014 Thermo Fisher Scientific Inc. All rights reserved. SilFlow is a registered trademark of SGE. ISO is a trademark of the International Standards Organization. All other trademarks are the property of Thermo Fisher Scientific Inc. and its subsidiaries. Specifications, terms and pricing are subject to change. Not all products are available in all countries. Please consult your local sales representative for details.

Africa-Other

+27 11 570 1840
Australia +61 3 9757 4300
Austria +43 1 333 50 34 0
Belgium +32 53 73 42 41
Canada +1 800 530 8447
China +86 10 8419 3588
Denmark +45 70 23 62 60

Europe-Other

+43 1 333 50 34 0
Finland/Norway/Sweden
 +46 8 556 468 00
France +33 1 60 92 48 00
Germany +49 6103 408 1014
India +91 22 6742 9434
Italy +39 02 950 591

Japan

+81 45 453 9100
Latin America +1 561 688 8700
Middle East +43 1 333 50 34 0
Netherlands +31 76 579 55 55
New Zealand +64 9 980 6700
Russia/CIS +43 1 333 50 34 0
South Africa +27 11 570 1840

Spain

+34 914 845 965
Switzerland +41 61 716 77 00
UK +44 1442 233555
USA +1 800 532 4752



Thermo Fisher Scientific,
 Austin, TX USA is
 ISO 9001:2008 Certified.

Thermo
 SCIENTIFIC

Part of Thermo Fisher Scientific